



CENTRIFUGES CS I / CS II

FLEXIBLE USE OF THE SYSTEM INCREASES LABORATORY EFFICIENCY INDIVIDUALLY CUSTOMIZABLE

CSI/CSII

Cytocentrifuges

CS I and **CS II** centrifuges are a cost-effective alternative for monolayer cell preparation, providing consistent performance and user friendliness. The systems are equipped with an open or closed rotor for up to 12 loading positions which can be easily removed for external loading.

Single or double sample chambers are available as disposable or reusable after proper cleaning. For a flexible use of the system a variety of sediment rotors are available on request. The **CS II** centrifuge is depending on the model equipped with cooling. Up to 10 memory programmes can be safed.

Highlights

Flexible selection and handling of rotors

- Cytocentrifugation open or closed (12 loading positions)
- Easily removable rotor for outside loading
- Large variety of rotors available (sediment, cytocentrifuge)
- CS II with optional cooling and up to 10 memory programmes

Suitable consumables

- Cytofunnel single or double (reusable or disposable)
- Large volume ECOfunnel® (reusable or disposable)
- Filter cards

Easy and safe handling

- Convenient preparation
- Slide clips applicable to reusable and disposable funnels
- · Slide label readable at any time
- Indication of the RCF during centrifugation

	Item No.		Item No.
CS I 230 V, 50/60 Hz Open rotor	15500000	CS II 230 V, 50/60 Hz Open rotor	15500100
CS I 230 V, 50/60 Hz Closed rotor	15500001	CS II 230 V, 50/60 Hz	15500101
CS I 115 V, 60 Hz	on demand	CS II 230 V, 50/60 Hz	
		Open rotor, with cooling CS II 230 V, 50/60 Hz	15501100
		Closed rotor, with cooling CS II 115 V, 60 Hz	15501101 on demand





CS rotor with 12 loading positions

	CSI	CSI	CSII	CSII
Standard components	Open rotor	Closed rotor	Open rotor	Closed rotor
Cellclip rotor,				
open (12 positions)	~		~	
Cellclip rotor,				
closed (12 positions)		~		•
Cellclip				
(stainless steel, 12 pcs.)	•	~	V	~
Cytofunnel,				
reusable (12 pcs.)	~		V	
Filter cards for cytofunnel,				
reusable				
Item No. 39910001 standard (200 pcs.)	v	~	~	•
Cellslides				
Item No. 39910003 (100 pcs.)	•	~	V	•
Mains cable	•	v	v	•
Manual	•	~	~	✓





Specifications	CSI	CSII
Application	cytology	cytology, laboratory
Cooling / temperature	-	down to -20 °C (optional)
Cytofunnel capacity	single 0.5 ml / chamber double 0.5 ml / chamber ECOfunnel® 5 ml / chamber	single 0.5 ml / chamber double 0.5 ml / chamber ECOfunnel® 5 ml / chamber
Memory programmes	-	up to 10
Rotor capacity	up to 12	up to 12
Applicable rotors	cellclip rotor (cytocentrifugation) swing-out rotor (different sizes) fixed angle rotor (different sizes)	cellclip rotor (cytocentrifugation) swing-out rotor (different sizes) fixed angle rotor (different sizes)
Centrifuge operation	up to 6,000 rpm depending on the selected rotor), timer 0 - 99 min.	up to 15,000 rpm depending on the selected rotor), timer 0 - 99 min.
Safety	automatic rotor recognition, lid locking and holding	automatic rotor recognition, lid locking and holding
Electrical parameters	230 V, 50/60 Hz	230 V, 50/60 Hz
Dimensions (W \times D \times H)	370 x 430 x 260 mm	400 x 700 x 350 mm
Weight	25 kg	29 kg


Reusable preparation systems

Cytofunnel single, reusable	Quantity	Item No.
Cellclip	1 pcs.	39910031
Cytofunnel single	12 pcs.	39910000
Filter cards (only for cytofunnel single 39910000) up to 0.5 ml	200 pcs.	39910001
Filter cards (only for cytofunnel single 39910000) thin, up to 0.4 ml	200 pcs.	39910006
Cellslides single, uncoated	100 pcs.	39910002
Cellslides single, coated	100 pcs.	39910003




Cytofunnel single, reusable

Cytofunnel double, reusable	Quantity	Item No.
Cellclip		39910031
Cytofunnel double	12 pcs.	
Filter cards for cytofunnel double up to 2 x 0.5 ml	500 pcs.	
Cellslides double, uncoated (only for cytofunnel double, reusable 39910010)	100 pcs.	39910012

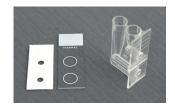

Cytofunnel double, reusable

Cytofunnel ECOfunnel®, reusable	Quantity	Item No.
Cellclip		39910031
Cytofunnel ECOfunnel®	12 pcs.	
ECOseal seal for ECOfunnel®	100 pcs.	
ECOslides, coated	100 pcs.	39910022

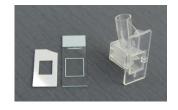

Cytofunnel ECOfunnel®, reusable

Accessories and consumables for third party products

All components of the CS centrifuge series like funnels, filter cards and slides are compatible to the Thermo Fisher / Shandon Cytospin series. To use CS centrifuge consumables in these third-party products, simply replace the original clip with the CS series cellclip Item No. 39910030.

Rotors

Rotor (closed)	1110.
CS centrifuge, removable, 12 positions 3991	0100



Rotor (open)	Item No.
CS centrifuge, removable, 12 positions	39910050



EASY rotor (closed)	item No.
10 marting with a trial and the claim	
12 positions, without stainless steel clips,	
for EASY / double cytofunnel	15500600



EASY rotor (open)	Item No.
12 positions, without stainless steel clips,	
for EASY / double cytofunnel	15500601



Swing-out rotor 8-point	Item No.
for 15 ml	39910042



Swing-out rotor 6-point	Item No.
for 50 ml	39910043
Adapter for 15 ml (only for 39910043)	39910044



The **cellclip of the CS series** an be easily exchanged for a third-party cellclip!



Cellclip of the CS series

Third-party cellclip

Disposable preparation systems

EASY cytology funnel can only be used in Thermo Cytospin or with the EASY rotor of the **CS** centrifuge!

EASY cytofunnel single, disposable	Quantity	Item No.
With white filter card up to 0.5 ml	40 pcs.	39910090
With brown filter card up to 0.5 ml		on demand





EASY cytofunnel single

EASY cytofunnel double, disposable	Quantity	Item No.
With white filter card up to 2 x 0.5 ml		39910091
With brown filter card up to $2 \times 0.5 \text{ ml}$		on demand





EASY cytofunnel double

Cytofunnel single, disposable	Quantity	Item No.
Cellclip	1 pcs.	39910031
Cytofunnel single up to 0.5 ml	100 pcs.	39910004
Cytofunnel single up to 0.5 ml		39910005
Cellslides single, uncoated	100 pcs.	
Cellslides single, coated	100 pcs.	39910003



Cytofunnel single, disposable

Cytofunnel double, disposable	Quantity	Item No.
Cellclip	1 pcs.	39910031
Cytofunnel double, sample chamber with 2 shifted holes, up to 2 x 0.5 ml	100 pcs.	39910014
Cytofunnel double, sample chamber with 2 shifted holes, up to 2 x 0.5 ml	500 pcs.	39910015
Cellslides double, uncoated (only for cytofunnel double 39910015)	100 pcs.	39910016
Cellslides double, coated (only for cytofunnel double 39910015)		on demand



Cytofunnel double, disposable

ECOfunnel®, disposable	Quantity	Item No.
Cellclip	1 pcs.	
ECOfunnel®, up to 5 ml	100 pcs.	
ECOfunnel®, up to 5 ml	500 pcs.	
ECOslides, coated		39910022
ECOslides, uncoated		on demand



ECOfunnel®, disposable

Service

We are committed to provide world-class support to our global customers at all times. We strive to advise you professionally before purchase and to support you during installation and initial setup. In addition, it goes without saying that we will always be available as a reliable partner during the longtime use of our products.

You can contact us directly with any kind of questions. In order to process your inquiries in time, please always include the following information:

- User name
- Address
- Phone number
- E-mail
- Type of device
- Serial number of device
- Date of purchase
- Detailed case description

Service Hotline: +49 (0) 6136 76997-0

Service E-mail: service@slee.de

Trainings

Service and application training

Well-trained staff is crucial for the implementation of regular service and maintenance of high quality.

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
We offer regular service and application training for our worldwide distribution partners.

Visit www.slee.de to access the training schedule and to download the registration form.

Training requests

Tel.: +49 (0) 6136 76997-0

E-mail: mail@slee.de





SLEE medical GmbH • Am Neuberg 14 55268 Nieder-Olm • Germany www.slee.de

T: +49 (0) 6136 76997-0 E: mail@slee.de









www.slee.de

YouTube

LinkedIn

Instagram

